

Lab 20 Discharge Device Upgrade

If a customer buys a new Lab 20 with the larger neck opening of 2.2 inches, they can keep their old manual discharge device provided they upgrade the rubber stopper. Please use part number 13717369 KIT DISCHARGE DEVICE UPGRADE. This kit includes a washer, larger rubber plug, and bolt assembly.



TEC 3000 – Changing the Logging Period

The default logging interval, rate at which data is recorded in the memory, of the TEC 3000 is 4 hours. Adjusting the logging period requires connecting the controller to the computer using a TEC COMKIT, Chart Connect 3000, and the ASCII command LOGPER. The logging period cannot be changed in the menu screens of the controller*. The TEC COM Kit can be ordered using part number PN: 13376947 and the Chart Connect program is available for download free of charge on the Chart Biomed website (<http://chartbiomed.com/tec2000.cfm>)

*NOTE: Changing the print interval only affects the controller's output when in printing mode.

TEC 3000 - Power Failure Alarm

Firmware versions before 2.00, displayed a power failure alarm if the battery voltage drops below 16 volts. Firmware version 2.00 and later added an additional parameter; a power failure alarm will be displayed if the battery voltage drops below 16 volts controller OR the controller has been operating using battery backup for 30 minutes.



8]gIf]Vi hYX`Vm



PrincetonCryo.com
Cryogenic Delivery and Storage Systems

www.PrincetonCryo.com | Sales@PrincetonCryo.com | 800.232.2796